

Hematoxylin and Eosin stain protocol for COT20

Time	Steps
120s	XTRA-SOLV
120s	XTRA-SOLV
120s	XTRA-SOLV
120s	EtOH 100%
120s	EtOH 100%
120s	EtOH 70%
120s	EtOH 50%
120s	H2O distilled
120s	HE Papanicolaou J.T.BAKER: 3873
120s	H2O tap
120s	*HCl Alcohol 0.2%
120s	H2O tap
120s	**H2O tap + NH4
120s	H2O distilled
120s	H2O distilled
120s	***Eosin 1% + Erythrosin B 1%
120s	***Eosin 1% + Erythrosin B 1%
120s	H2O tap
120s	EtOH 96%
120s	EtOH 100%
120s	XTRA-SOLV
120s	XTRA-SOLV
52min	